

Genorise Red Blood Cell Lysis Solution

Genorise Red Blood Cell Lysis Solution is to isolate white blood cells from 0.3 ml of blood for isolations of DNA or protein, and is for 1000 applications. If a smaller or larger sample volume is employed, the reagent quantity should be proportionally decreased or increased. This kit will guarantee the quality and quantity of white blood cells and is much cheaper than the similar products (\$68/1000 ml).

Materials provided in the kit:

1000 ml Red Blood Cell (RBC) Lysis Solution

Materials required but not provided:

PBS, pH 7.4, prepared in sterile distilled deionized water.

Protocol

Red Blood Cell Lysis

1. Take 0.3 ml whole blood to a 1.7 ml microcentrifuge tube, and add 0.9 ml of Red Blood Cell Lysis Solution, vortex for 20 sec.
2. Incubate at room temperature for 10 min to completely disrupt the red blood cells.
3. Centrifuge the tube at 5000 x g for 5 min at 4°C and discard the supernatant.
4. Add 1 ml of cold PBS to the white pellet, suspend the cells by a pipette, and centrifuge the tube at 5000 x g for 5 min at 4°C.
5. If the white pellet contains red substance, add 0.3 ml RBC Lysis Solution, suspend the cells by a pipette, and incubate at room temperature for 5 min.
6. Centrifuge the tube at 5000 x g for 5 min at 4°C and discard the supernatant.
7. Repeat Step 5 and 6 until the pellet is red-free.